


Since 1954, SPEX CertiPrep is the industry leader in the CRM marketplace meeting the needs of laboratories worldwide with innovation and research. Accredited by A2LA to ISO/IEC 17025:2017 & ISO 17034:2016. Certified by DQS to ISO 9001:2015.

Part #: ODER-250G  
Size: 250 g

Directions for use: Open container and place in affected area (lid can be removed by cutting off). As the pellets are exposed to air contaminants, the purple pellets will turn brown. Shake material occasionally to expose more of the purple pellets. When a majority of the pellets have turned brown, it is time to replace the product. Depending on exposure, the product will last up to 2-3 months. To reorder, call 1-800-LAB-SPEX or 732-549-7144 or visit [www.spexcertiprep.com](http://www.spexcertiprep.com)

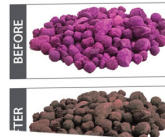


## OdorEroder

Neutralizes Odor-Causing Chemicals

SPEX CertiPrep®  
203 Norcross Avenue, Metuchen, NJ 08840  
732-549-7144 • [www.spexcertiprep.com](http://www.spexcertiprep.com)

Part #: ODER-100G  
Size: 100 g

Directions for use: Open container and place in affected area (lid can be removed by cutting off). As the pellets are exposed to air contaminants, the purple pellets will turn brown. Shake material occasionally to expose more of the purple pellets. When a majority of the pellets have turned brown, it is time to replace the product. Depending on exposure, the product will last up to 2-3 months. To reorder, call 1-800-LAB-SPEX or 732-549-7144 or visit [www.spexcertiprep.com](http://www.spexcertiprep.com) (Part #ODER-100G).

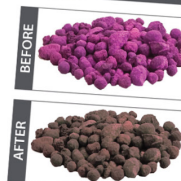

## OdorEroder®

### Contamination Control Products

OdorEroder® effectively neutralizes a wide range of offensive chemical odors and fumes in the lab, everything from Aldehydes to Xylenes. These non-toxic, environmentally safe purple beads are placed where odor-causing chemicals in the air pass near the beads. When this occurs, the purple beads absorb and chemically transform the odor-causing chemicals into harmless compounds that remain trapped within the beads. As the purple beads absorb, they start turning brown. When a majority of the beads have turned brown, it is time to replace your OdorEroder®. Depending on exposure, the OdorEroder® lasts up to three months. Contact [CRMSales@spex.com](mailto:CRMSales@spex.com) for your FREE sample. For additional product information and for a FREE sample of the OdorEroder®, please visit [www.spexcertiprep.com](http://www.spexcertiprep.com).

### OdorEroder® is Effective in the Following Areas:

- Hoods
- Lab Benches
- Glove Boxes
- Lab Refrigerators
- Waste Disposal Areas
- Chemical Storage Cabinets
- Other Odor-Causing Areas Within the Lab

Advantages of OdorEroder® Versus Carbon	OdorEroder®	Carbon
Highly effective at neutralizing volatile compounds	√	—
Absorbs and chemically transforms chemical odors into harmless compounds that remain trapped within the product	√	—
Changes color to signal replacement of the product	√	—

Ordering Information	
Description	Part Number
OdorEroder - 100 g	ODER-100G
OdorEroder - 250 g	ODER-250G

### Contact Us

Phone: 800.LAB.SPEX • 732.549.7144 • Fax: 732.603.9647  
[CRMSales@spex.com](mailto:CRMSales@spex.com) • [spexcertiprep.com](http://spexcertiprep.com)

© 2020 SPEX CertiPrep. All Rights Reserved.

CONNECT WITH US

